



## PuraLev® Life Science Pump Series

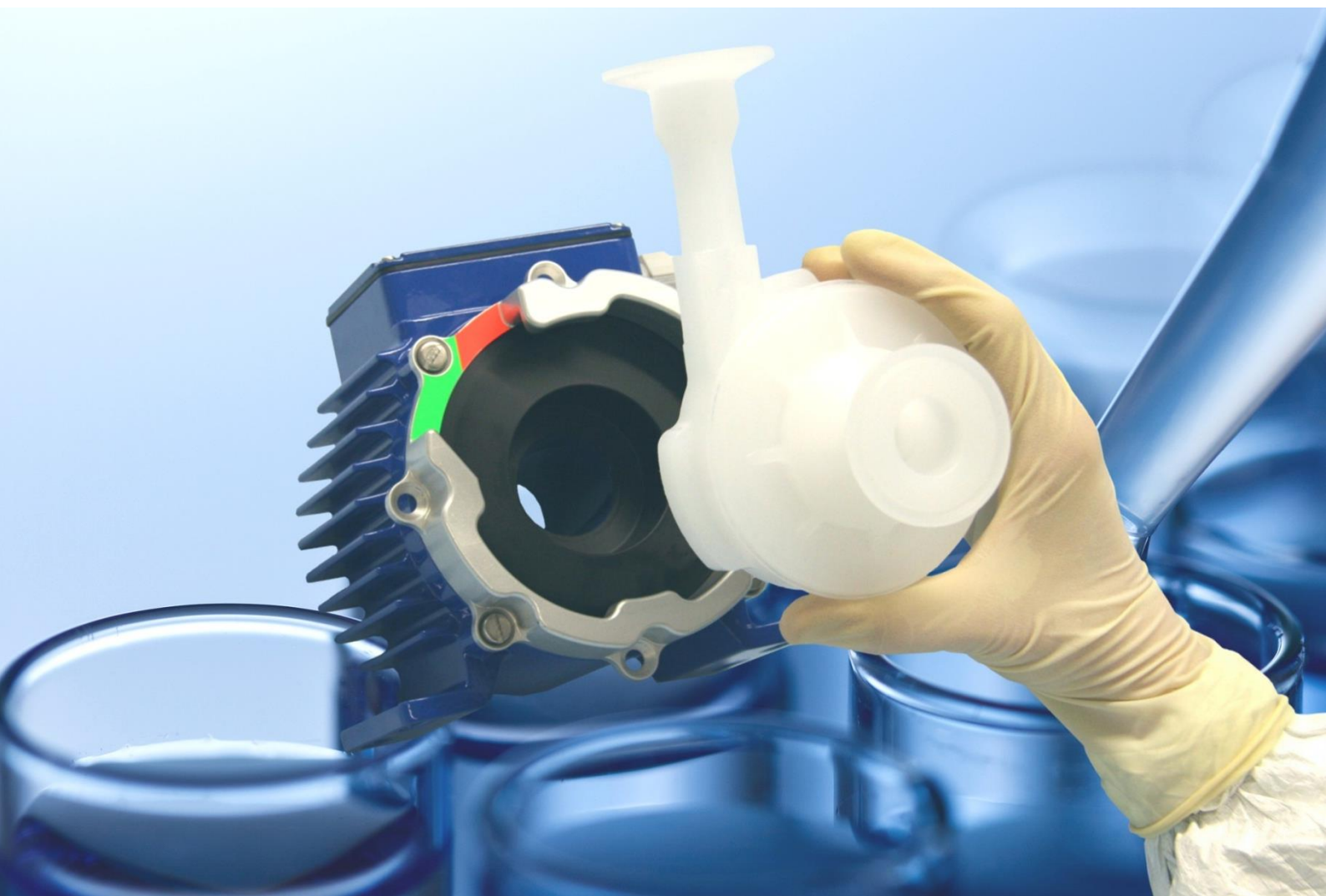


### PuraLev® 600SU (Single-Use)

3.1 bar	(45 psi)
75 liters/min	(20 gallons/min)


Low Shear Design - High Cell Viability

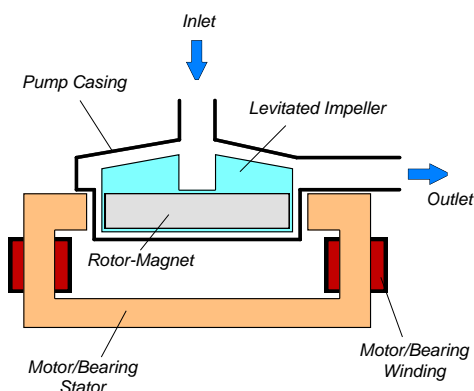

**Figure 1:** Schematic of the main elements of the maglev centrifugal pump

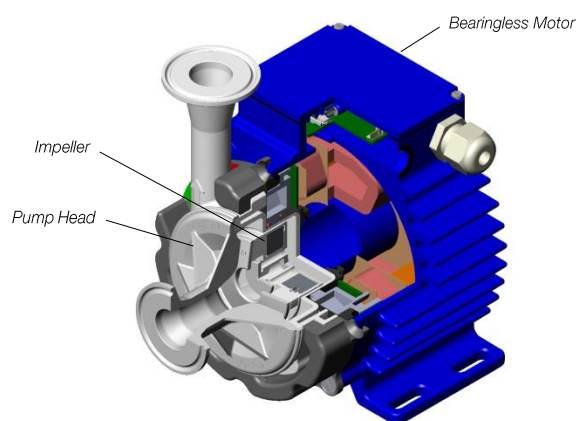

**Figure 2:** Cross-section of the bearingless pump motor and pump head

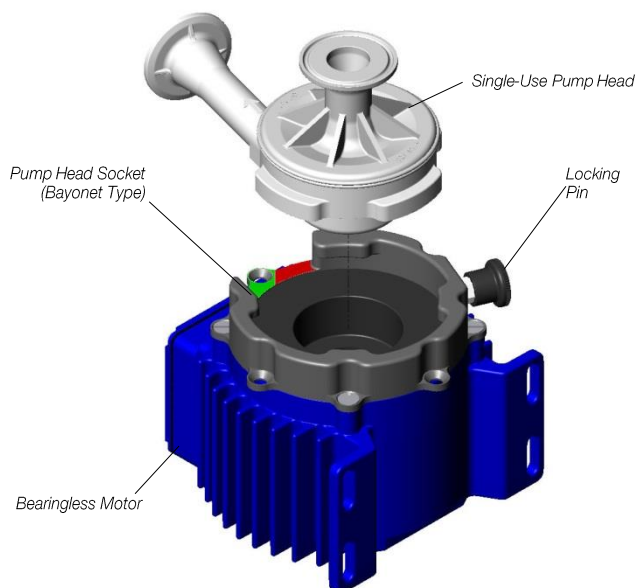

**Figure 3:** Single-use pump head concept

## INTRODUCTION

Levitronix® has developed a revolutionary pump that has no bearings to wear out or seals to break. Based on the principles of magnetic levitation, the pump's impeller is suspended, contact-free, inside a sealed casing and is driven by the magnetic field of the motor (Figure 1). The impeller and casing are both fabricated from biocompatible (FDA, USP-VI, BSE/TSE and Animal free) gamma sterilizable polypropylene (PP) and together they make up the disposable pump head. A simple and intuitive exchange of the single use pump head is achieved with a bayonet socket type mounting procedure (see Figure 5). Flow rate or pressure are precisely controlled by electronically regulating the rotor speed, which eliminates any pulsation. With the lack of mechanical bearings plus the self-contained pump head design, the risk of contamination is drastically reduced. The absence of narrow gaps between the impeller and pump casing, plus the low-shear pump design allows the gentle pumping of sensitive liquids. The pump casing is fabricated with Triclamp fittings and can be easily inserted and removed with an intuitive bayonet socket.

## SYSTEM BENEFITS

- Low shear-forces
- Reduced risk of contamination due to the self-contained design with magnetic bearings
- No particle generation
- No over-pressure situations (compared to roller pumps)
- No narrow gaps between the impeller and pump casing where bacteria could be entrapped
- Pump head is gamma sterilizable
- Biocompatibility of wet materials: FDA, USP-VI, Animal/BSE/TSE free
- Bayonet socket design for easy and intuitive exchange of disposable pump head (see Figure 5)
- Small size
- Dry running capability
- Proven technology in the medical (disposable blood pumps) and semiconductor (high-purity pumps) industries
- High flow capability with compact design
- Pulsation free

## APPLICATIONS

- Pumping of shear-sensitive liquids and cells
- Bioprocessing (for example perfusion)
- Recirculation and transfer applications in bioreactors
- Filtration

## STAND-ALONE SYSTEM CONFIGURATION

The stand-alone configuration of the *PuraLev® 600SU* pump system consists of a controller with an integrated user panel allowing the operator to set the speed manually (see *Figure 6*). The speed is automatically stored in the internal EEPROM of the controller. As an option, the speed can also be set with an analogue signal (see specification for *Position 3a* in *Table 2*).

## EXTENDED SYSTEM CONFIGURATION

The extended version of the *PuraLev® 600SU* pump system (*Figure 7*) consists of a controller with an extended PLC interface. The PLC interface allows the speed to be set via an external signal, facilitating precise closed-loop flow or pressure control when either a flow or pressure sensor is integrated into the system (see specification of *Position 3b* in *Table 2*). A computer can be connected via a USB interface to allow communication with *Levitronix® Service Software*. Hence parameterization, firmware updates and failure analysis are possible.

## ATEX / IECEx SYSTEM CONFIGURATION

An ATEX / IECEx certified motor together with the pump head allows installation of motor and pump head within an ATEX Zone 2 area (see *Figure 8*). The ATEX / IECEx motor (*Pos. 2b* in *Table 2*) comes with special connectors and relevant extension cables (*Pos. 5a* and *5b* in *Table 3*). An Ex conform solution is needed for the motor cables to leave the ATEX area. One option is an ATEX certified cable sealing system as listed in *Table 4* (see *Pos. 9*) and shown in *Figure 12*.

- ATEX / IECEx certified for Category 3G and 3D (Zone 2 for Gas and Zone 22 for Dust).
- Thermal classification T4 (< 110 °C = 230 °F) for maximum liquid temperature of 90 °C / 194 °F.
- ATEX / IECEx marking of motor with pump head:
 

CE Ex II 3G Ex c nAc IIC 110°C (T4)  
 CE Ex II 3D Ex c tc IIIC T110°C IP67
- Explosion groups:
 

Group IIA: Propane (IPA), Methane, Acetone, Acetaldehyde  
 Group IIB: Ethylene, Ethylenglycol  
 Group IIC: Acetylene, Hydrogen (not carbon disulphide)
- ATEX / IECEx listing corresponds to UL hazardous location Class 1 Division 2.

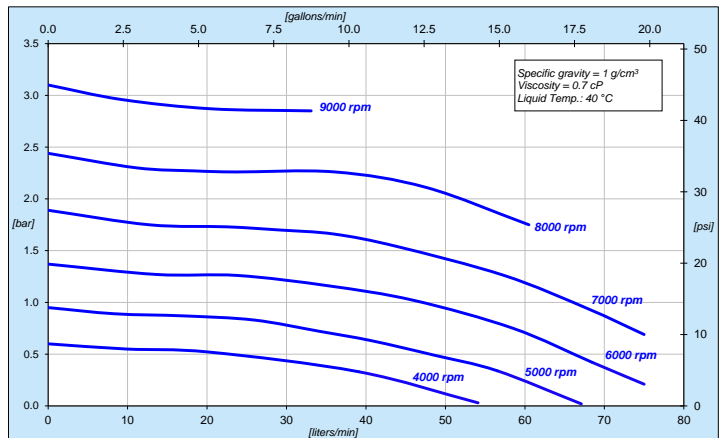


Figure 4: Pressure/flow curves (DCP-600.2 pump head)




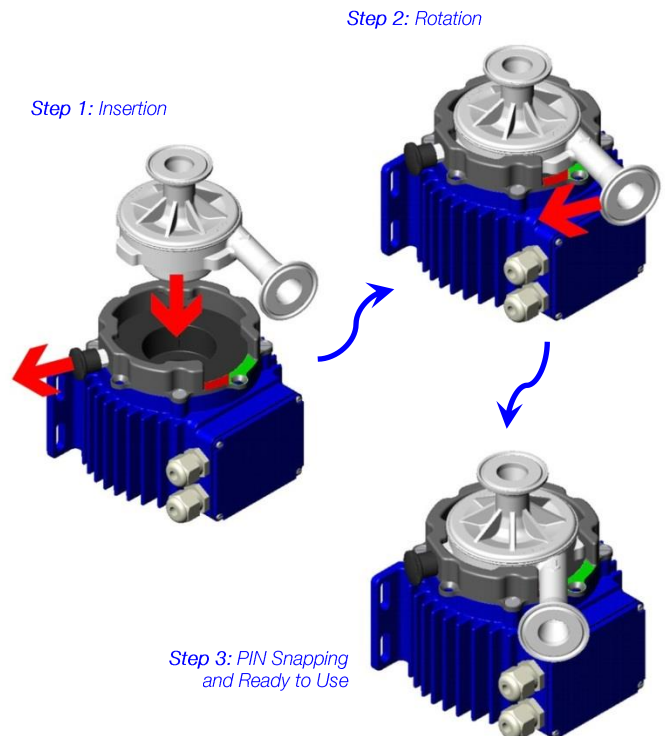



Figure 5: Intuitive 3-step pump head mounting procedure with bayonet type socket (PHS-600.1)

# SYSTEM CONFIGURATIONS

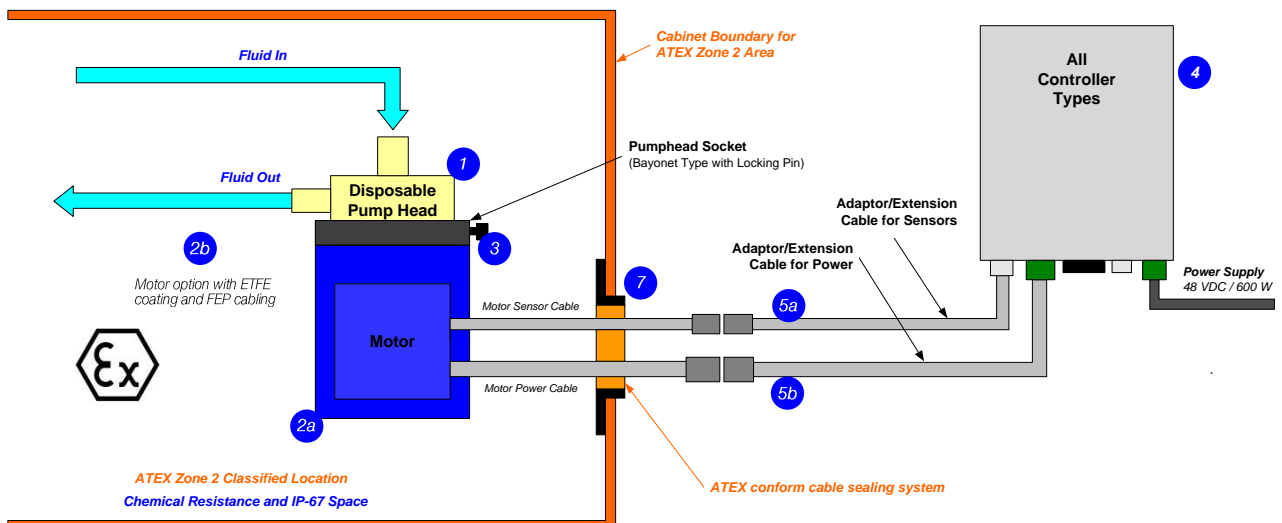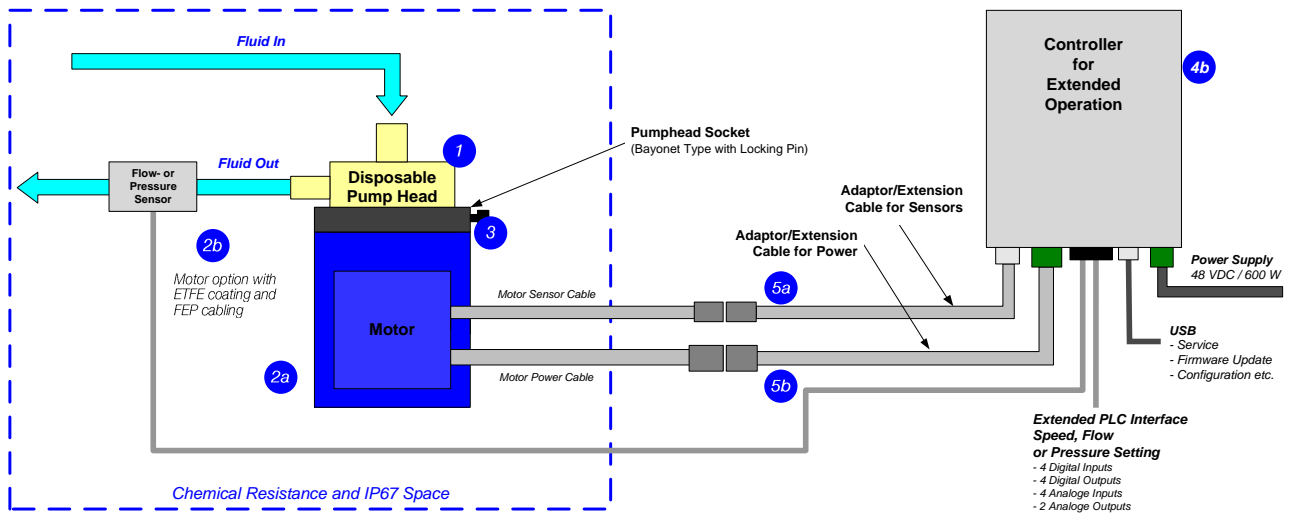

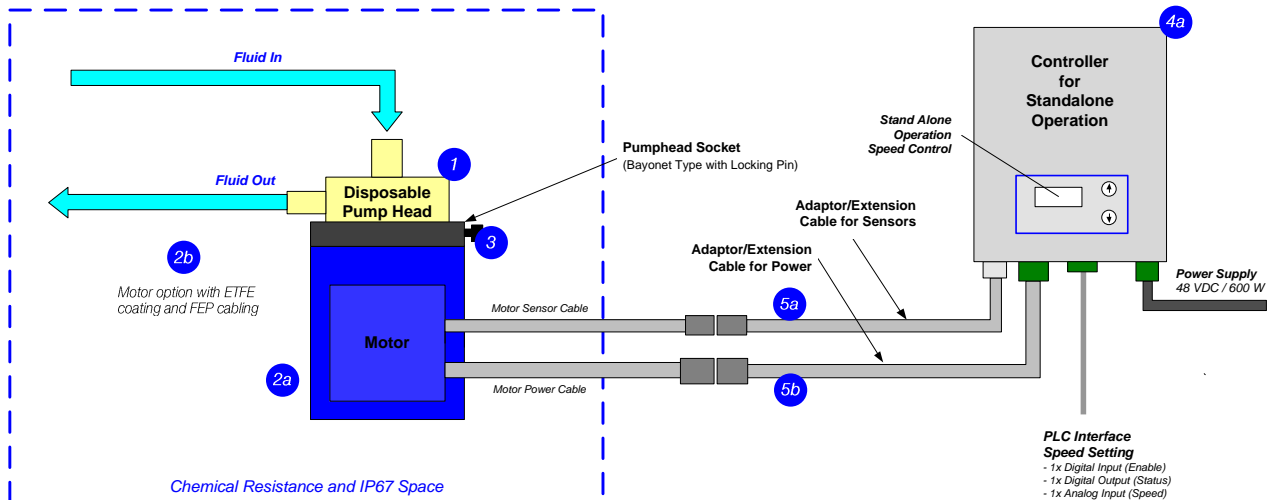

## DIMENSIONS OF MAIN COMPONENTS

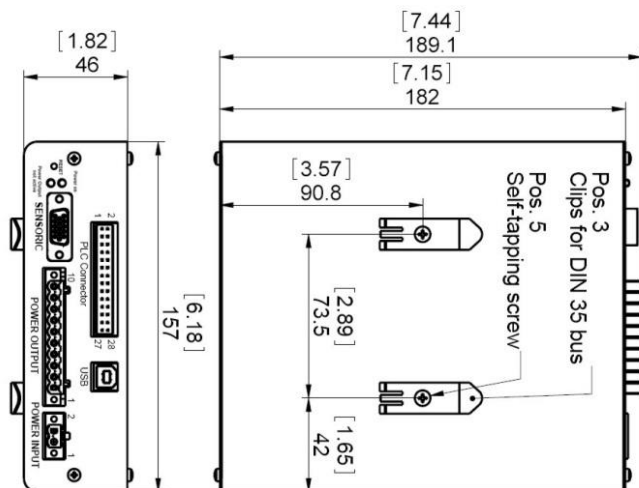

Controller  
LPC-600.1-01

4a





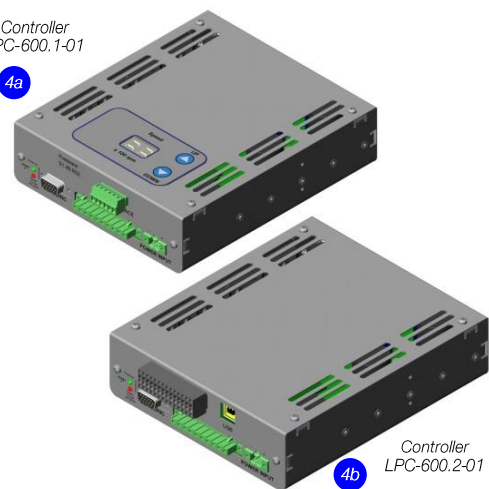


Figure 9: Dimensions of controllers

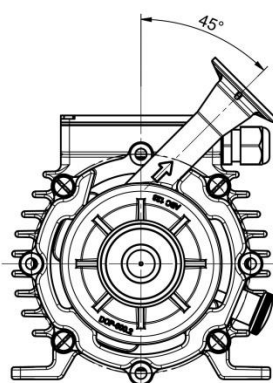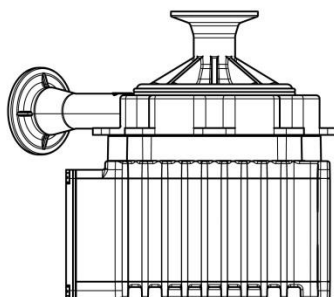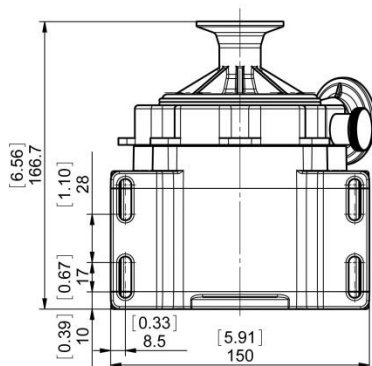

Standard Mounting Position

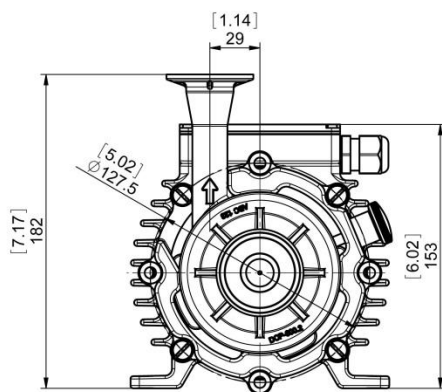

Vertical Mounting Position

Pump Head Socket  
PHS-600.1


3

Single Use Pump Head  
DCP-600.1

1

Inlet/Outlet Fitting:  
1" Triclamp  
ASME BPE 2009  
ID = 22.1mm

Snap-PIN for  
Pump Head Fixation

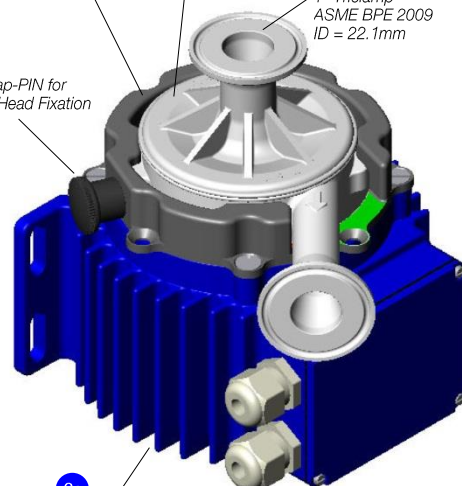

Motor  
LPM-600.5

2a

Figure 10: Dimensions of motor with single-use pump head

## ORDER INFORMATION

System Name	Article #	Pump Head Socket	Motor	Controller	Note
PLD-600SU.1	100-90654	PHS-600.1	LPM-600.5	LPC-600.1-01	Adaptor/Extension (0.5 - 10m) cables according to Table 3 have to be ordered as separate article with specified length. ATEX Cable Sealing System can be ordered according to Table 4. Certifications: CE, IECEx CB scheme, ETL (NRTL), ATEX and IECEx. <sup>1</sup>
PLD-600SU.2	100-90655		LPM-600.5	LPC-600.2-01	
PLD-600SU.4	100-90657		LPM-600.4	LPC-600.1-01	
PLD-600SU.5	100-90658		LPM-600.4	LPC-600.2-01	

**Table 1:** Standard system configurations with motor, pump head socket and controller  
1: Certified components are available on request.

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature
1a	Single-Use Pump Head	DCP-600.2	100-90784	Impeller / Pump Housing Housing Sealing In-/Outlet Fittings Max. Flow Max. Diff.-Pressure Max. Viscosity Wet Pump Volume/Surface Max. Liquid Temp. Sterilization Methods	Polypropylene (FDA, USP Class VI, BSE/TSE/Animal free) Infrared welding Triclamp 1" (ASME BPE 2009) 75 liters/min / 20 gallons/min 3.1 bar / 45 psi 50 cP 113 ml / 370 cm <sup>3</sup> 60°C / 140°F Gamma radiation up to 40 kGy.
1b	Irradiated Pump Head	DCP-600.2-G25	100-91079	Applied Gamma Dosage	≤25 kGy
2a	Motor (ATEX / IECEx)	LPM-600.5	100-10039	Housing Cable / Connectors ATEX / IECEx Marking	Epoxy (anti-corrosive) coated Aluminum, waterproofed (IP67) 2x 3m cables with PVC jacket / 2x circular (M23, IP-67) CE II 3G Ex c nAc IIC T110°C (T4) CE II 3D Ex c tc IIC T110°C IP67
2b	Motor (ATEX / IECEx)	LPM-600.4	100-10038	Housing Cable / Connectors	ETFE (chemical resistant) coated Aluminum, waterproofed (IP67) 2x 3m cables with FEP jacket / 2x circular (M23, IP-67)
3	Pump Head Socket	PHS-600.1	100-90696	Mounting Type Material Assembly Screws	Bayonet type with locking pin Anodized Aluminum 4 pcs M6x16 mm (Stainless Steel)
4a	Standalone Controller (User Panel)	LPC-600.1-01	100-30039 (Power supply cable and Enable connector incl.)	Voltage / Power Interfaces Standard Firmware	48V DC / 600 W Panel to set speed (automatic storage on internal EEPROM) PLC with 1x analog input ("Speed") 4 - 20 mA 1x digital input ("Enable") 0 - 24 V (optocoupler) 1x digital output ("Status") 0 - 24 V (relais) D9.25
4b	Extended Controller (PLC and USB)	LPC-600.2-01	100-30040 (Power supply cable and PLC connector incl.)	Interfaces Standard Firmware	PLC with - up to 4 digital inputs 0 - 24V (optocoupler) - up to 4 digital outputs 0 - 24 V (relais) - up to 2 analog inputs 4 - 20mA - up to 2 analog inputs 0 - 10 V - up to 2 analog outputs 0 - 5 V USB interface (for service and system monitoring) D9.48

**Table 2:** Specification of standard components

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature
5a	Extension Adaptor Cable for Sensors	MCAS-600.2-05 (0.5m) MCAS-600.2-30 (3m) MCAS-600.2-50 (5m) MCAS-600.2-70 (7m) MCAS-600.2-100 (10m)	190-10226 190-10238 190-10127 190-10105 190-10239	Jacket Material Connectors	PVC Circular Wallmountable, Metallic (IP-67) to D-SUB
5b	Extension Adaptor Cable for Power	MCAP-600.2-05 (0.5m) MCAP-600.2-30 (3m) MCAP-600.2-50 (5m) MCAP-600.2-70 (7m) MCAP-600.2-100 (10m)	190-10227 190-10240 190-10126 190-10106 190-10241	Jacket Material Connectors	PVC Circular Wallmountable, Metallic (IP-67) to COMBICON

**Table 3:** Specification of adaptor/extension cables

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature
6	Air Cooling Module	ACM-600.2	190-10140	Material / Connection Port Air Pressure	PP (+ 40% Talkum) / NPT 1/4" ~1 - 3 bar (14 - 43 psi)
7 (a - f)	ATEX Cable Sealing System	ACS-A.1 (Roxtec)	100-90292	Sleeve (a) and Gasket (b) Frame (c) 2x Cable Module (d)	Stainless Steel and EPDM Roxylon (EPDM rubber) Roxylon (EPDM rubber)
8	AC/DC Power Supply	TSP 600-148-M (M = Modified Levitronix design from Traco)	100-40013 (Traco ID Number: T1068-01A)	Voltage / Power Output Voltage Input Certification or Standards	48 VDC / 600 W 85 - 265 VAC (automatic detection) CB, UL, CSA, Semi F47

**Table 4:** Specification of accessories

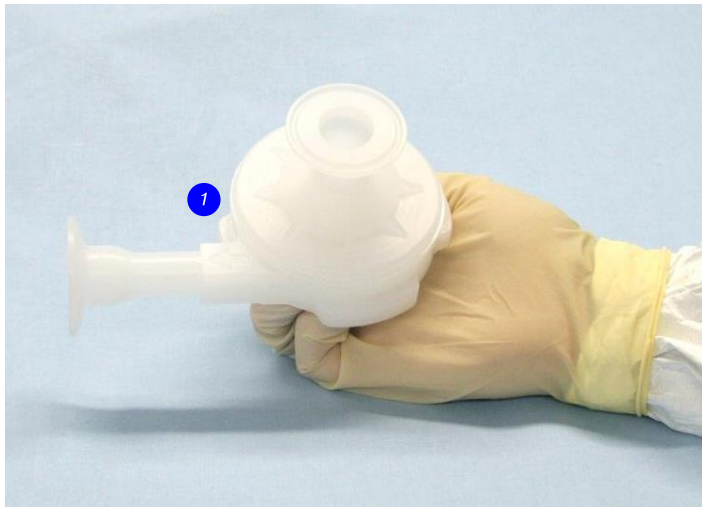
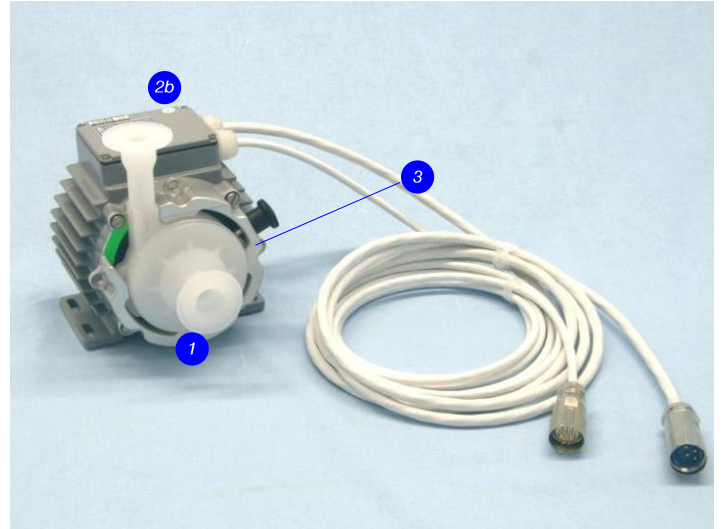


Figure 11: Pump system with standard components

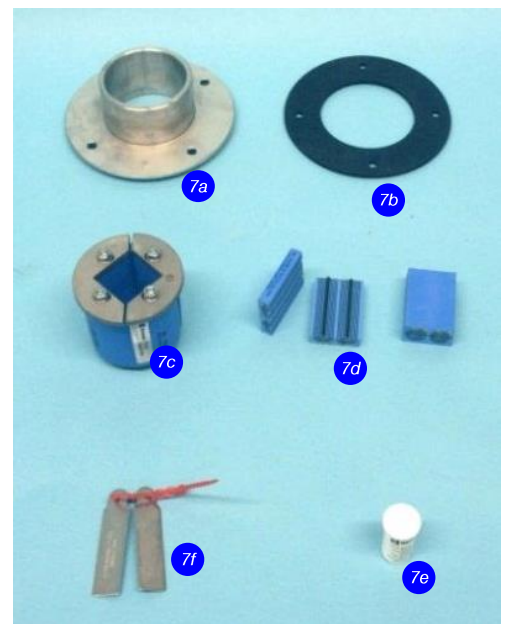


Figure 12: Accessories

*Levitronix®* is the world-wide leader in magnetically levitated bearingless motor technology. *Levitronix®* was the first company to introduce bearingless motor technology to the Semiconductor, Medical and Lifescience markets. The company is ISO 9001 certified. Production and quality control facilities are located in Switzerland. In addition, *Levitronix®* is committed to bring other highly innovative products like the *LEVIFLOW®* flowmeter series to the market.



#### Headquarter and European Contact

Levitronix GmbH  
Technoparkstr. 1  
CH-8005 Zurich  
Switzerland

Phone: +41 44 445 19 13  
Fax: +41 44 445 19 14  
E-Mail: [salesEurope@levitronix.com](mailto:salesEurope@levitronix.com)

#### US Contact

Levitronix Technologies LLC  
20 Speen Street, Suite 102  
Framingham, Massachusetts 01701  
USA

Phone: +1 508 861 3800  
Fax: +1 508 861 3837  
E-Mail: [salesUS@levitronix.com](mailto:salesUS@levitronix.com)

#### Japan Contact

Levitronix Japan K.K.  
Wing Eight 5floor, 4-16-4  
Asakusabashi, Taito-ku  
Tokyo, 111-0053 Japan

Phone: +81 3 5823 4193  
Fax: +81 3 5823 4195  
E-Mail: [salesJapan@levitronix.com](mailto:salesJapan@levitronix.com)

#### Taiwan Contact

Levitronix Taiwan  
5F, No. 251, Dong Sec. 1, Guangming  
6th Rd., Chu Pei City,  
Hsin-Chu 302, Taiwan, R.O.C.

Phone: +886 3 657 6209  
Fax: +886 988 321472  
E-Mail: [salesAsia@levitronix.com](mailto:salesAsia@levitronix.com)